
LAB TEACHER DEMO WORKSTATION

1114K, 1114KF, 1116K, & 1116KF SERIES ASSEMBLY INSTRUCTIONS

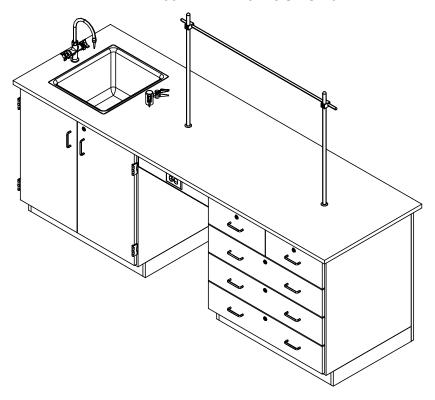

TOOLS REQUIRED
CONTACT CEMENT
SCREW DRIVER
C-CLAMPS
SHIM STOCK (IF NEEDED)

ASSEMBLY COMPONENTS					
ITEMS INCLUDED	PART#	PART DESCRIPTION	QTY		
TOP	VARIES	TOP, (STYLE VARIES), 1.00X30X96,BLK, (EITHER FIXTURE OR FLAT)	1		
CABINET	N/A	B86 - 36"W X 35"H X 26.25"D	1		
CABINET	N/A	B23 - 36"W X 35"H X 26.25"D	1		
BACK PANEL	N/A	94"W X 35"H X 0.75"D	1		
APRON	N/A	0.81"THK X 4.50"W X 22"L	1		
SCREWS - BACK PANEL	100478	SCREW,#8X1 ¼ SQ. WASHER FLAT HEAD	12		
L BRACKET	100355	BRACKET, RAIL 1.25X1.25X4.25	2		
SCREWS - L BRACKET	100649	SCREW,#8X¾ PHIL TRUSS HEAD SMS	8		
SILICONE	100711	SILICONE,SEALANT,CLEAR	1		
RUBBER BASE MOLDING	100283	RUBBER BASE MOLDING (SOLD BY FOOT)	25		
STAINLESS STEEL CORNERS	100057	CORNER-STAINLESS STEEL, BASE	8		
SCREWS - BASE	100064	SCREW,#6X½ PTH SMS SS SCREW	32		
ELECTRICAL BOX	100033	ELECTRICAL BOX 3.00X2X2.5 DEEP	1		
ELECTRICAL OUTLET	100034	ELEC,OUTLET,GFI,20A 125V, BLK	1		
FIXTURE - FAUCET	100048	FIXTURE,L4100FH-132AWSA 90 DEGREE-G	1		
FIXTURE - FAUCET	100467	FIXTURE,L414 HOT & COLD MIXING	1		
FIXTURE - SINK	100699	SINK,EPOXY L30,BLK	1		
FIXTURE - SINK TRAP	100056	SINK TRAP,PLASTIC ADJUSTABLE	1		
FIXTURE - SINK STOPPER	100112	SINK STOPPER 1 ½,BLK	1		
FIXTURE - STRAINER	100055	SINK,OUTLET/STRAINER EPOXY,BLK	1		

NOTE:

- 1. TOPS ARE EITHER PHENOLIC OR EPOXY.
- 2. **XX** INDICATES ITEMS USED ONLY FOR THE "K" SERIES.




ASSEMBLY COMPONENTS				
ITEMS INCLUDED	PART#	PART DESCRIPTION	QTY	
ROD SOCKET	206505	ROD SOCKET ASSY,SOCKET,WASHER,NUT	2	
CROSSBAR	100006	rod,crossbar ¾ dia.x48 l	1	
UPRIGHT	100004	ROD,UPRIGHT ¾ DIA.X36 L	2	
CLAMP	100002	ROD CLAMP ¾ X ¾ ASSY	2	

NOTE:

- 1. TOPS ARE EITHER PHENOLIC OR EPOXY.
- 2. **XX** INDICATES ITEMS USED ONLY FOR THE "K" SERIES.




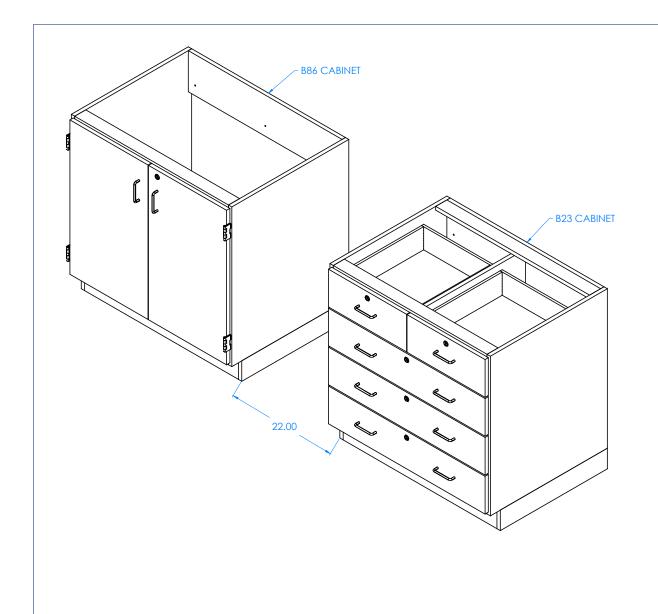

STEP 1:

- POSITION THE TWO CABINETS IN DESIRED LOCATION.
- IF YOU HAVE A "K" SERIES UNIT, PLACE UNIT OVER UTILITY HOOKUPS.
- LEAVE ENOUGH ROOM IN BETWEEN CABINETS FOR APRON.
- LEVEL CABINETS.
- ATTACH TO FLOOR.




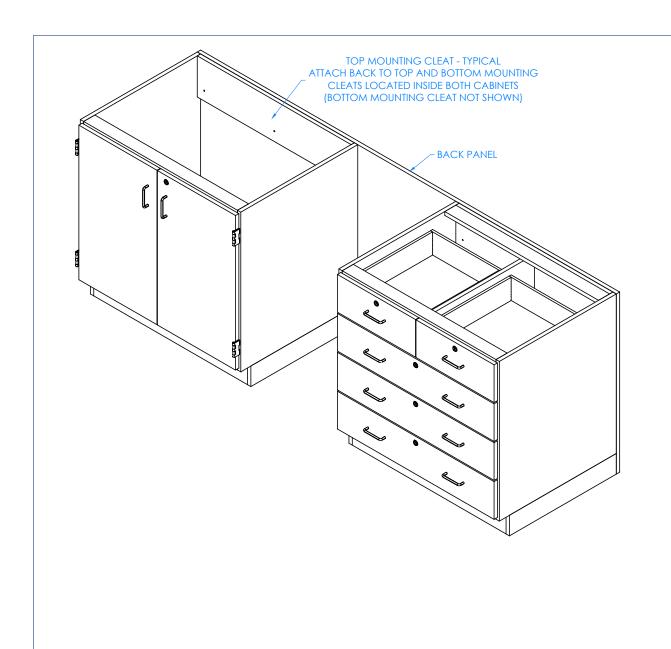

STEP 2:

- ATTACH BACK PANEL TO BACK OF CABINETS WITH HARDWARE PROVIDED.
- SCREW INTO TOP AND BOTTOM MOUNTING CLEATS LOCATED INSIDE BOTH CABINETS.
- MAKE SURE BACK PANEL IS FLUSH WITH OUTSIDE EDGES OF CABINETS.




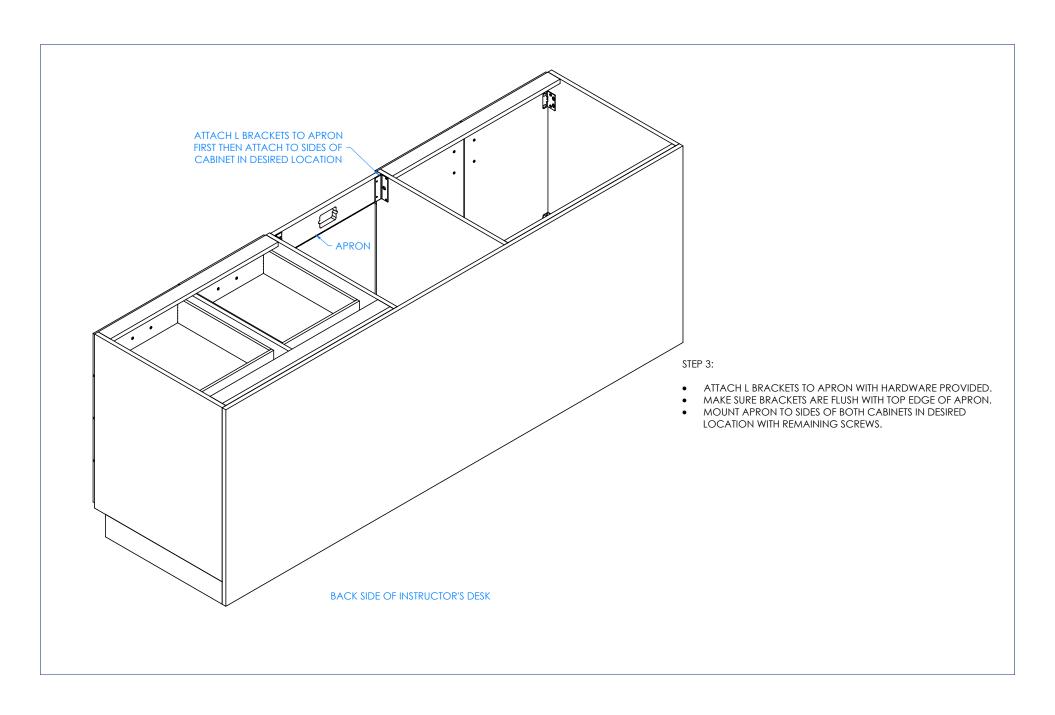



STEP 4:

- APPLY SILICONE ADHESIVE TO TOP EDGES OF CABINETS AND APRON.
- PLACE TOP ON UNIT WITH PROPER OVERHANG (1" OVERHANG ON SIDES AND 1.50" ON FRONT AND BACK).
- CLAMP ON TOP TO TOP RAILS OF CABINETS AND APRON TO REMOVE ANY WARP USING SHIMS AND C-CLAMPS, IF NEEDED, WHEN ALIGNING THE TOP (DO THIS BEFORE ADHESIVE ON EDGES SETS UP).
- LET CURE FOR 6-12 HOURS.

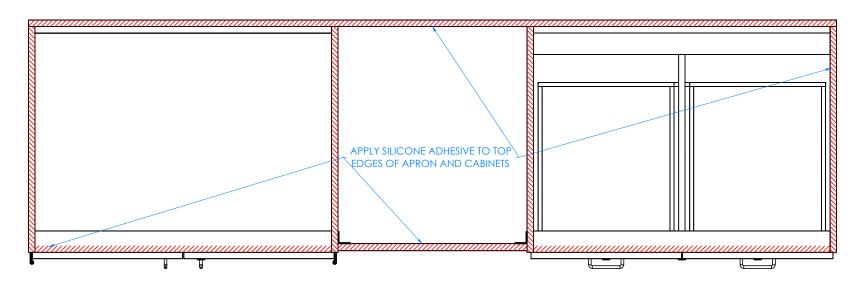


TOP VIEW OF INSTRUCTOR'S DESK SHOWN




STEP 5:

- IF YOU DO NOT HAVE A "K" SERIES UNIT, SKIP TO STEP 6.
- APPLY SILICONE ADHESIVE TO SINK OPENING ON TOP.
- LOWER SINK INTO OPENING.
- WIPE OFF ANY EXCESS CAULK AND ALLOW TO SET.
- LET CURE FOR 6-12 HOURS.

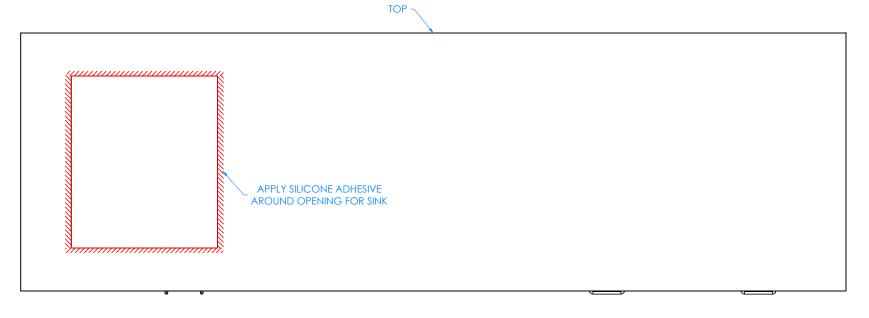

TOP VIEW OF INSTRUCTOR'S DESK SHOWN



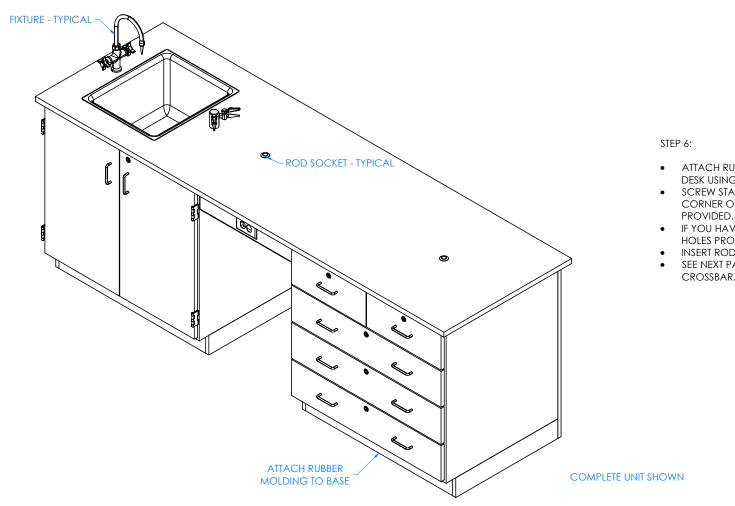


- ATTACH RUBBER MOLDING TO BASE OF INSTRUCTOR'S DESK USING CONTACT CEMENT.
- SCREW STAINLESS STEEL CORNER BRACKETS ON EACH CORNER OVER BASE MOLDING WITH HARDWARE PROVIDED.
- IF YOU HAVE A "K" SERIES UNIT, INSTALL FIXTURES IN THE HOLES PROVIDED ON TOP.
- INSERT ROD SOCKETS INTO PRE-DRILLED HOLES IN TOP.
- SEE NEXT PAGE ON HOW TO INSTALL UPRIGHTS AND CROSSBAR.


NOTE:

 ALL WATER, ELECTRICAL, AND GAS CONNECTIONS SHOULD BE PERFORMED BY A TRAINED PROFESSIONAL PER LOCAL CODE.




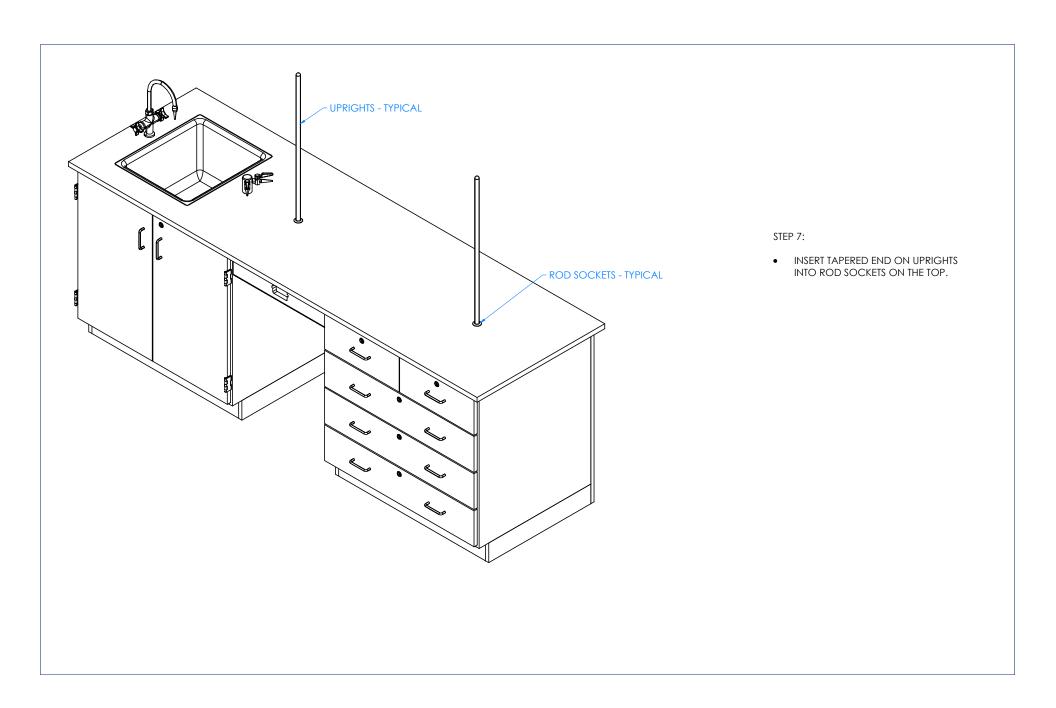




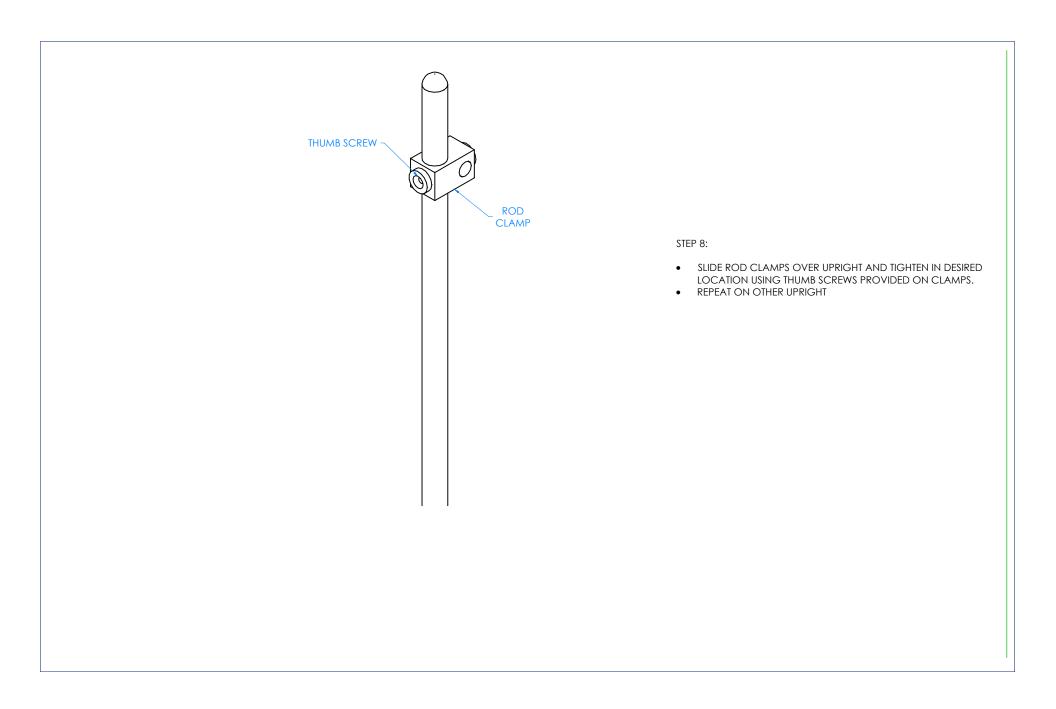



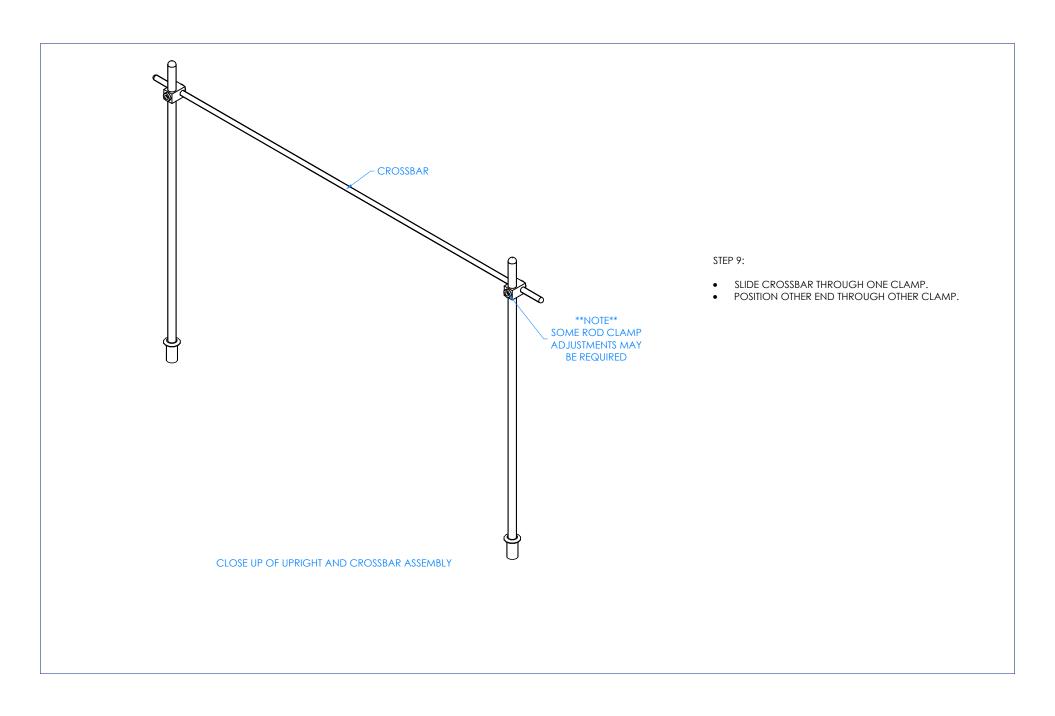



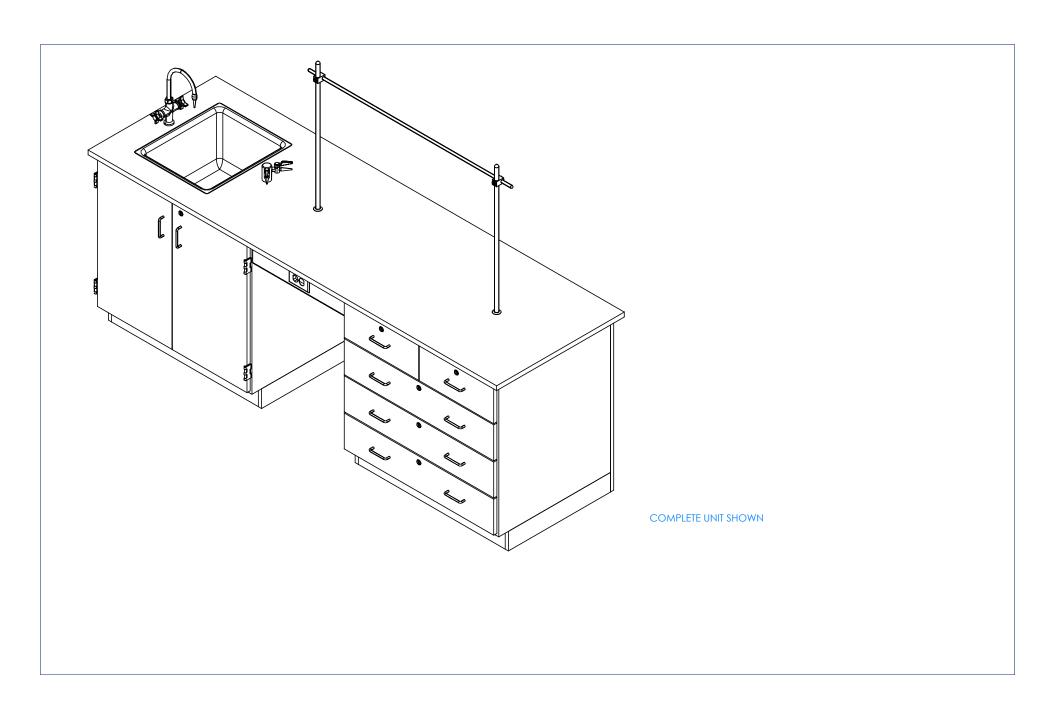



TROUBLESHOOTING YOUR GFCI ELECTRICAL RECEPTACLE:

<u>PLEASE NOTE: THROUGHOUT ANY OF THE FOLLOWING STEPS (PROCEDURES); IF YOU ARE NOT SURE YOU CAN DO THIS JOB SAFELY, AND COMPETENTLY, REFER THIS WORK TO QUALIFIED PROFESSIONAL!</u>

If your unit comes equipped with a GFCI (Ground Fault Circuit Interrupter) electrical receptacle, and there is no power in the receptacle please consider/check the following:

- The extension cord (if being used) is working properly.
- The breaker hasn't been tripped, or turned off.

This type of GFCI has two testing-related buttons on it. One button is appropriately labeled "TEST", and the other button is labeled "RESET". To test the GFCI receptacle follow these steps:

- Plug in an appliance (lamp or night light) into the outlet. The light should now be on. Then press the "TEST" button on the GFCI. The GFCI "RESET" button should pop out, and the light should go out.
- If the "RESET" button pops out, but the light doesn't go out, the GFCI has been improperly wired. In this case please contact a certified professional. There may also be a problem with other wiring in the same circuit.
- If the "RESET" button doesn't pop out, the GFCI is defective, or malfunctioned, and should be replaced.
- If the GFCI is functioning properly, and the lamp goes out, press the "RESET" button to restore power to the outlet.

Conversely, if you have a GFI that has tripped (which is common) and it will not reset, you may have a wiring short in the circuit, a defective appliance on the circuit, or the GFI itself has become defective. To test a tripping GFCI follow these steps:

- Remove every appliance connected to the GFCI's circuit and reset it. If it doesn't reset there may either be a wiring fault behind a socket outlet, or your GFCI itself has become faulty.
- Make sure whatever you are plugging into to the GFCI is dry and not damaged.
- Only plug in one item at a time. If you are plugging in a defective item it will cause the GFCI to trip, and that item therefore should be replaced.

If you are still having difficulty the easiest way to troubleshoot a GFI is to obtain a GFI tester, which is available at most hardware stores. It plugs into the GFI outlet, and will supply you with a "snapshot" of your connections, indicating wiring problems and/or the condition of the GFI. Another way to troubleshoot is to simply purchase a new GFI and install it.

